

# LOUISIANA DEPARTMENT OF TRANSPORTATION & DEVELOPMENT VERIFICATION PROCEDURE

## WATER BATH

### EQUIPMENT

1. Ruler
2. Secondary Calibrated Thermometer

### PROCEDURE

1. Measure and record the distance from the perforated shelf to the bottom of the water bath.
2. Measure and record the distance from the bottom of the bath to the surface of the water.
3. Measure and record the temperature of the water in the bath. Temperature may be +/- 0.1°C of the set point.

EQUIPMENT VERIFICATION RECORD

Inspected By: \_\_\_\_\_ Date \_\_\_\_\_

Apparatus: **WATER BATH (T245, TR 305)** Make: \_\_\_\_\_

Identification No.: \_\_\_\_\_ Verification Freq. 12 MO.

Previous Verification Date: \_\_\_\_\_ Next Due Date: \_\_\_\_\_

Verification Equipment Used: \_\_\_\_\_

Verification Procedure: \_\_\_\_\_

Distance from perforated shelf to bottom of bath (min. 2"): \_\_\_\_\_

Distance from bottom to surface of water (min. 5½"): \_\_\_\_\_

Bath thermostatically controlled? (yes or no): \_\_\_\_\_

Temperature of water (140 ± 1°F.: 60 ± 0.5°C.): \_\_\_\_\_

Action Recommended: Repair \_\_\_ Replace \_\_\_ None \_\_\_

Comments \_\_\_\_\_

CC: ENGINEERING TECHNICIAN MANAGER, EQUIPMENT TESTING